DOOR ALARM

Por favor lea las instrucciones atentamente antes de usar!
Please read the instructions carefully before use!

Precauciones/Caution

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Pharmacy Refrigerator

ENGLISH

This **Operation Manual** applies to the following models: TH-076 | TH-081 | TH-176

Table of Contents

1. Application Notes	26
2. Safety Instructions	27
3. Precautions in Use	29
4. Product Installation	29
4.1 Installation Enviroment	29
4.2 Installation Site.....	29
4.3 Preparation before use.....	29
4.4 First Power-on	30
4.5 Operation after Power Failure.....	30
5. Product composition and overview	32
6. Operating Instructions	34
6.1 Function Introduction	34
6.2 Optional function.....	41
7. Maintenance and Service	41
7.1 Equipment Maintenance	42
7.2 Equipment Discontinuation	42
7.3 Maintenance, Replacement and Recovery of Rechargeable Batteries	42
8. Troubleshooting and Maintenance Services	43
9. Specifications	44
10. Packaging List	44

1. Application Notes

Thank you for choosing and using Thermolabil® products! For your safety and convenient use and reasonable maintenance of the equipment, please read the Operating Instructions carefully before use and keep them properly for reference.

The equipment operator can copy some chapters of this operation manual, but only for internal use, for example, to instruct the user how to deal with emergencies. These chapters are clearly marked in the catalogue of the manual.

Thermolabil® has no obligation or responsibility for instrument damage caused by the user's failure to use the equipment according to the instructions or method specified by the manufacturer.

Due to the rapid improvement of Thermolabil® products, the functions described in the instructions may be inconsistent with those of the products you purchased. Please refer to the physical functions.

- Please read carefully the Attention and Safety Precautions (in 2. Safety Instructions).
- During transportation or use, no violent vibrations or collisions are allowed and the refrigerator shall be kept away from rain. Store in a clean room with humidity no more than 80 %, no corrosive gas and good ventilation.
- Pharmacy/medical refrigerators (hereinafter referred to as equipment) can only be operated by trained and authorised personnel.
- Maintenance of the equipment can only be completed out by Thermolabil® or an agent authorized by Thermolabil®.
- If the operator encounters any situation not mentioned in this manual, please contact Thermolabil® or the agent authorised by Thermolabil® for the correct handling method.
- If the equipment is not used according to the method specified in the specifications, it may be damaged.
- Try to use the accessories provided by Thermolabil®. If users would like to use other accessories, Thermolabil® will not be responsible for the adverse consequences caused therefrom.
- Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.

Tips

- ✓ Properly use protective equipment (including protective clothing, protective gloves, goggles, etc.).
- ✓ Keep good hygiene habits.
- ✓ Please pay attention to safety when using this product.
- ✓ Keep clear of obstruction all ventilation openings in the appliance enclosure or in the structure for building-in.
- ✓ Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.
- ✓ Do not damage the refrigerating circuit.
- ✓ Do not use electrical appliances inside the food/ice storage compartments unless they are of the type.
- ✓ Recommended by the manufacturer.

2. Safety Instructions

When using this product for the first time, please pay attention to the meaning of the following warning signs and carefully read the safety precautions, so that you can use this equipment safely and correctly.

 Warning! Failure to observe precautions may result in serious personal injury or death.	 Attention! Failure to observe precautions may result in personal injury or equipment failure and related material loss.
 Protective conductor terminal.	 Risk of Explosion This signal indicates the risk of explosion when using volatile and explosive chemicals.
 Beware of fire. Warning; flammable material	 Warning: crushing of hands

Manufacturer

Series NO

Date of production

 Warning: Failure to observe the precautions may result in serious personal injury or death.
 Do not touch the equipment with wet hands to avoid electric shock; Equipment shall not be grounded through gas pipeline, water supply pipeline, telephone line or lightning rod, which may easily cause electric shock.
 This equipment can only be installed by professional technicians or after-sales service personnel, it may cause electric shock or fire.
 Be sure to install the equipment on a solid and flat ground and take due care to prevent tipping over. If the ground is not solid enough or the installation location is not appropriate, it may cause the equipment fall over and cause damage or personal injury.
 Please handle the power cord carefully to avoid short circuits or open circuits. Please turn off the power pulling out the power plug. Hold the power plug carefully and pull it out. Do not pull the wires of the power plug. Otherwise, it may cause electric shock or fire due to short circuit. Do not bundle the power cord, do not press it under furniture or heavy objects, and do not get close to heat sources such as compressors.
 Please insert the power plug into the outlet tightly to ensure firmness and reliability, so as to avoid electric leakage; After installation, the power plug must be within reach, so that the power plug can be unplugged in time in case of emergency.
 Separate special outlets must be used and grounded reliably. The cross-sectional area of the copper conductor in the wall connected with the outlet must be more than 4 mm ² . Do not lengthen the power cord without authorisation to avoid heating or fire.
 Do not use the power supply that is not specified in the equipment design, so as to avoid overheating, short circuit and other faults. For example, connecting 110 V rated voltage products to 220 V power supply may cause faults such as overheating and equipment burning. For detailed input voltage AC, please refer Specification (rated voltage ±10%). If the operating voltage is too low or too high, a suitable automic voltage stabiliser must be installed for cooperative use.
 Please place the equipment stably and avoid shaking.
 Do not place the equipment in a dangerous area, and do not operate the equipment near flammable items to prevent explosion or fire accidents.

	Do not place the equipment in areas exposed to sun or rain, so as to prevent danger such as short circuits or overheating.
	Do not tilt or lay the equipment sideways, and do not impact the equipment body ; Refrigeration systems are installed in the equipment, which is easy to damage by tilt or impact.
	Please place the equipment in a dry and dust-free environment to avoid risks such as overheating and short circuit.
	In case of unexpected sound, smell, smoke, etc...when power is turned on, please unplug the power in time and contact the manufacturer or supplier.
	Please place the equipment in a dry and ventilated environment, and ensure that the equipment vents and instrument surfaces are not blocked or shielded by walls or other objects. Do not use it in a poorly ventilated environment to prevent damage caused by the heat released by equipment.
	It is forbidden to disassemble and modify this equipment without authorisation, so as to avoid possible safety hazards. In this case, Thermolabil® will not bear any responsibility for quality accidents.
	Está prohibido colocar las mercancías peligrosas inflamables y explosivas, ácidos corrosivos fuertes, álcalis en el equipo y otros artículos que son inadecuados para el equipo.
	It is forbidden to put inflammable and explosive dangerous goods, strong corrosive acids, alkalis and other items that are unsuitable for the equipment in the equipment.
	Metal objects such as nails or iron wires shall not be inserted into any aperture and gap or any outlet of the equipment, otherwise electric shock or injury may be caused to accidental contact between the above objects and moving parts.
	In order to ensure the normal operation and ventilation and heat dissipation of the equipment, left and right sides of the cabinet shall be at least 30 cm away from the wall, and the air inlet and air outlet must not be blocked by obstacles.
	This equipment must be connected to a ground wire

	Note: Failure to observe the precautions may result in personal injury or equipment failure and related property losses.
	It is forbidden to store living animals, flowers or other items with strict temperature requirements in the equipment.
	When the equipment is running, do not touch the inner surface of the cabinet without wearing protective gear.
	Hold the handle and close the door to avoid pinching your fingers; When the equipment is not in use for a long time, please unplug it and pack it for storage.
	When restarting the equipment after a power failure or power off, please check the equipment settings, otherwise, the stored items may be damaged due to the change of settings.
	The equipment can be used for item preservation, not as production equipment.
	Keep the keys properly, so as to avoid accidents when children accidentally open the door accidentally.
	When handling the equipment, please be careful not to tip it over to the equipment, so as to prevent equipment damage or personal injury.
	When handling, it shall be lifted from the bottom, with the inclined plane not be greater than 45°, and it shall be handled with care. Please use the equipment in safe areas. Improper use harm to human health or the environment.

3. Precautions in Use

- Before putting the items into the equipment, please confirm that the temperature in the freezer storage chamber has reached the set value and then put the articles in batches. Every time you put in items, they shall not exceed 1/3 of the inner volume of the freezer storage chamber so as to prevent excessive temperature rise.
- The equipment temperature display value is the temperature at the temperature sensor in the freezer storage chamber. There is a certain gap between the displayed temperature and the actual temperature at the centre of the equipment when the equipment just starts running, but as the equipment enters a stable state, the displayed temperature will gradually approach the actual temperature.
- Please use a diluted neutral cleaning to clean the equipment, and do not use brushes, acid, gasoline, soap powder, polishing agent or hot water to clean the equipment, otherwise the painted surface and plastic rubber parts may be damaged. Be careful not to wipe plastic rubber parts with volatile solvents such as gasoline.

4. Product Installation

4.1 Installation Environment

- + Ambient temperature: 16°C ~ 32°C, the most ideal temperature is 18°C ~ 25°C, and the air conditioning system shall be used when necessary.
- + Relative humidity: ≤ 80%RH.
- + There is no strong vibration and corrosive gas around.
- + Avoid the existence of a large amount of dust.
- + Avoid rocking or shaking the equipment.
- + Elevation of the working position of the equipment: less than 2000m.
- + Indoor use, pollution degree 2 and overvoltage category II.
- + For detailed input voltage AC, please refer to Specification (rated voltage ±10%).
- + There is no direct sunlight, other cold and heat sources and strong electromagnetic interference, which will affect the normal operation of the control system and directly damage the system in severe cases.

4.2 Installation Site

In order to operate the equipment normally and obtain the best performance level, the installation site of the equipment shall meet the following requirements:

- + It cannot be installed in a narrow and closed space, and the door of the room shall not be smaller or lower than this equipment, so as to avoid maintenance difficulties in case of equipment failure in case of equipment failure, which may result in damage to stored items due to the failure to repair the equipment in time;
- + The installation floor must be solid, flat, non-combustible and able to bear the weight of the equipment during operation;
- + It shall be with good ventilation, and direct sunlight shall be avoided;
- + Each equipment needs to use a power outlet independently. Please ensure that the plug and outlet are firmly connected;
- + Check the working voltage before use. In areas with unstable voltage, consider using a voltage stabiliser suitable for equipment load to ensure that the input voltage requirements in the installation environment are met;

Attention: Since the ambient temperature has a great influence on the equipment, if the above environmental requirements cannot be met, the equipment may not run normally. Please improve the environment before using the equipment; the equipment is operated intermittently.

4.3 Preparation before use


- 1 Remove the outer packaging of all products (including the protective foam in the packaging box).
- ⚠ **Warning:** Don't put the plastic bags within the reach of children, so as to prevent suffocation accidents.
- 2 Inventory of accessories: Please check the accessories and materials according to the packing list.

- 3 Cleaning: Clean the product once before use.
- 4 Before use, remove the small wire shelf and put it on the bottom of the cabinet (applicable to the following models: TH-076).
- 5 Installation of door handle. In order to install handles, the handles shall be placed together with accessories.

Handle installation method:

Firstly, open the door handle as shown in the following figure; Then, remove the screw for installing the door handle on the cabinet and install the handle seat with this screw; Finally, install the handle cover, insert both ends of the handle cover into the corresponding grooves of the handle seat and then press it to close it.

- 6 Installing the fixing plates
 - a Firstly take out the two fixing plates and check if consistent.
 - b Remove the bolts from the fixing plate, and then fix the plate on the back of the refrigerator.
 - c Conect the fixing plate to the immovable wall or bracket on the back of the refrigerator and finally check whether it is fixed.

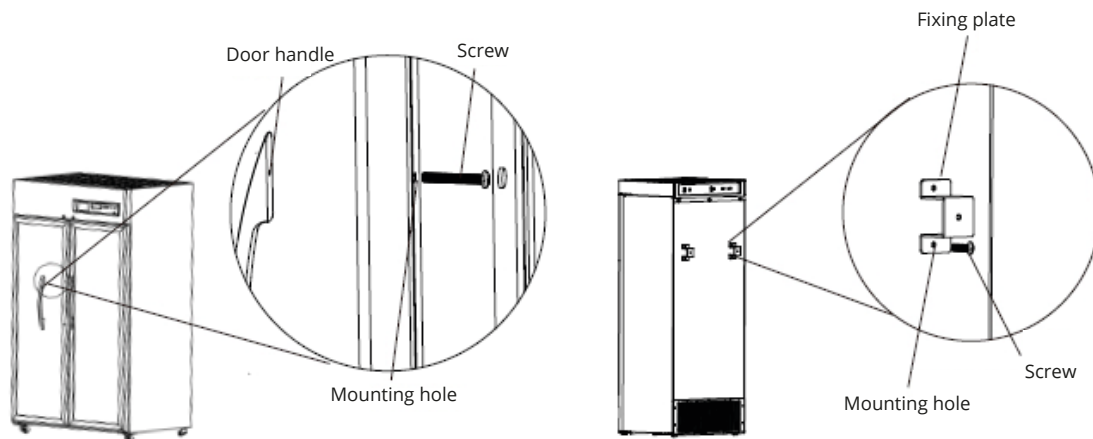4.4 First Power-on

When using the equipment for the first time, please follow these steps:

- 1 After the equipment is placed, leveled and cleaned, it shall stand more than 24 hours and then power it on to ensure the normal operation of the equipment.
- 2 Under no-load conditions, connect the power cord to a special outlet with the appropriate specifications.
- 3 After powering on, turn on the power switch of the equipment. Toogle the on/off switch to the ON position located on the back of the unit.
- 4 Check whether the operating temperature of the equipment reaches the required value, observe the normal start and stop of the equipment for more than 24 hours, and put a small amount of items in the refrigerator after confirming the normal performance.
- 5 Please store items in batches, with items not exceeding 1/3 of the cabinet volume each time. Ensure that the equipment is runnig properly after shutdown for more than 12 hours before putting in the next batch of items).
- 6 Try not to open the door during cooling, otherwise the temperature will rise.

4.5 Operation after Power Failure

- 1 The equipment has a memory function for the set value. When the power is restored after power failure, the equipment will continue to operate before power failure.
- 2 Once the equipment is powered off, it requires 5 minutes before it can be turned on again, so as to avoid damaging the compressor.
- 3 We guarantee the normal operation of this equipment under certain conditions, but we are not responsible for any loss or damage of stored items after power failure.

! Notes:

- A special person shall be responsible for checking and recording the running status of the equipment every day (record and check once every 2-4 hours). In case of failure or shutdown, the temperature in the refrigerator will rise. If it cannot be repaired in a short time, please take out the stored items and transfer them to a place that meets the temperature requirements for storage to avoid damage to the items.
- Before putting items into the equipment, it shall be confirmed in advance whether the temperature range of the equipment meets the temperature requirements of the items, so as to avoid damage to stored items due to the difference between the settable temperature of the equipment and the required temperature of the items. Please pay attention not to block the air outlet and air inlet when putting items into the equipment.
- Due to refrigeration inertia, there is a certain difference between the actual display temperature and the set temperature of the equipment, which is a normal phenomenon.
The equipment is an item storage equipment, which cannot be used for routine production operations.
- It is strictly forbidden to put too many items which are relatively hot into the equipment at the one time, otherwise the compressor will run for a long time, and be burned due to high temperature. Items must be put in batches, so as to ensure that the refrigerator is cooled step by step until the temperature required for storing items is reached.
- Electrical appliances without production licence shall not be used inside the equipment.
- Do not change the set temperature frequently in a short time, otherwise, the expected setting effect may not be achieved due to the large temperature inertia; Ensure that there is certain air circulation space around the cabinet when putting items, especially do not block the temperature sensor in the cabinet (for collecting the temperature of the cabinet), otherwise it will affect the stability and accurate control of the temperature in the cabinet.
- Items shall not be placed directly at the bottom of the refrigerator, but on the bottom shelf, otherwise the refrigeration effect of the equipment will be affected.
- When putting in items, if the moisture content of the items is too much or too little, it will affect the humidity change in the cabinet, so it is best to keep the items sealed; The humidity of the working environment will affect the change of humidity in the cabinet, especially if the door is opened frequently and the door is not closed properly.

! Warning:

- Children are not allowed to play with this equipment as a game prop, otherwise the injury or loss caused therefrom will be at their own risk.
- To avoid the risk of electric shock, this equipment must only be connected to a mains supply with protective earth.
- No modification of this equipment is allowed.
- Do not modify this equipment without authorization of the manufacturer.
- Protection impairment if used in a manner not specified by the manufacturer.
- The instructions concerning persons (including children) with reduced physical, sensory or mental capabilities, or lack of experience and knowledge, or children playing with the appliance are not required.
- Keep all ventilation openings in the enclosure or, in the structure for building in structure, clear of obstructions.
- Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.
- Do not damage the refrigerant circuit.
- Protection impairment if used in a manner not specified by the manufacturer.

5. Product composition and overview

Image 1. Refrigerator TH-076

*The **Refrigerator TH-076** contains three shelves, each with a capacity of 13kg.

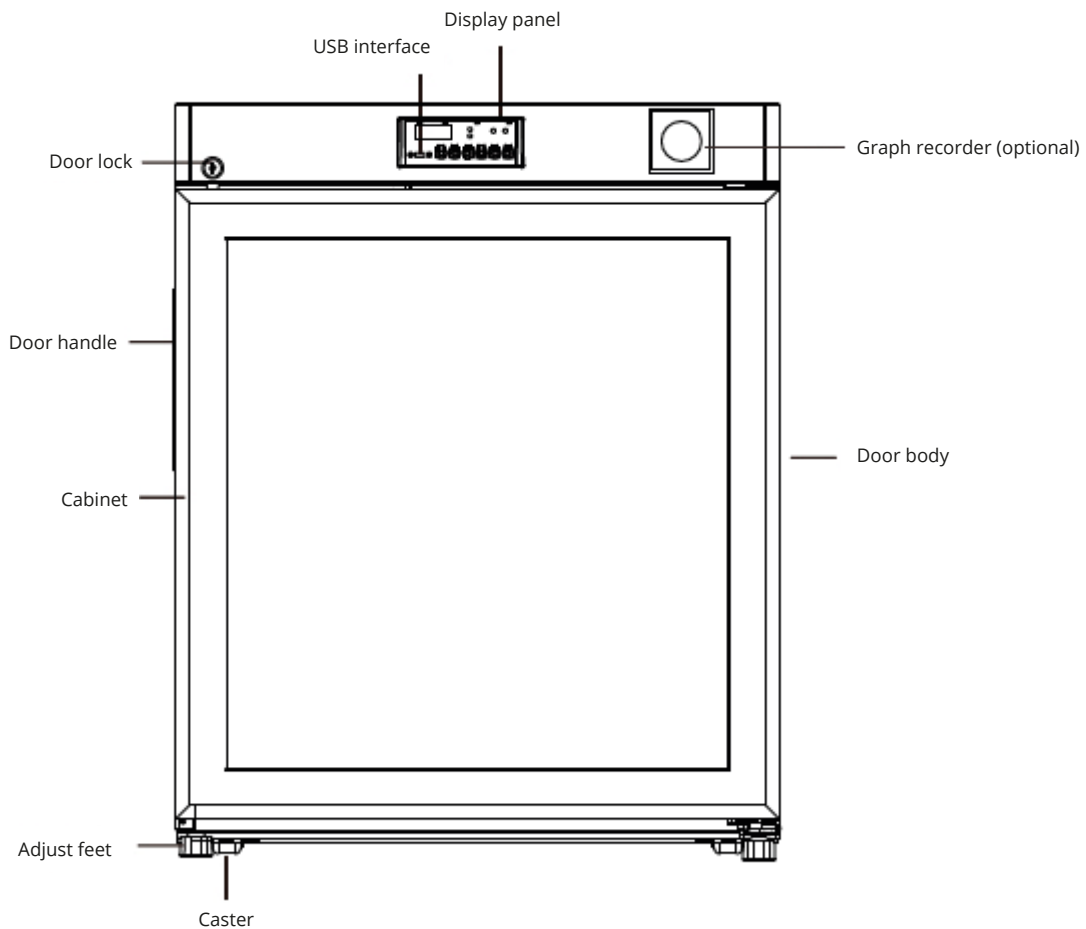
Image 2. Refrigerator TH-081

*The **Refrigerator TH-081** contains two shelves and a suspension frame, each with a capacity of 21Kg.


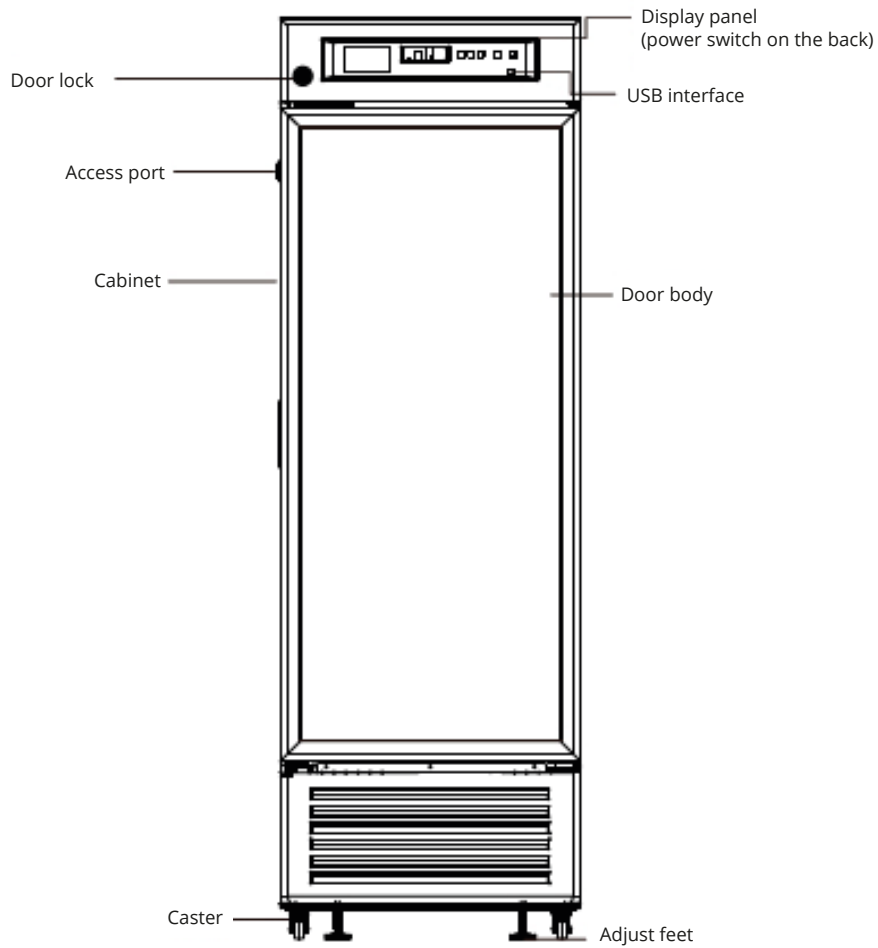
Image 3. Refrigerator TH-176

*The **Refrigerator TH-176** contains four shelves and one wire basket, each with a capacity of 25 kg.

6. Operating Instructions


6.1 Function Introduction

This series include two control systems, based to which different adjustment methods apply. Please select the proper adjustment method according to the type of control system with the storage box you purchased.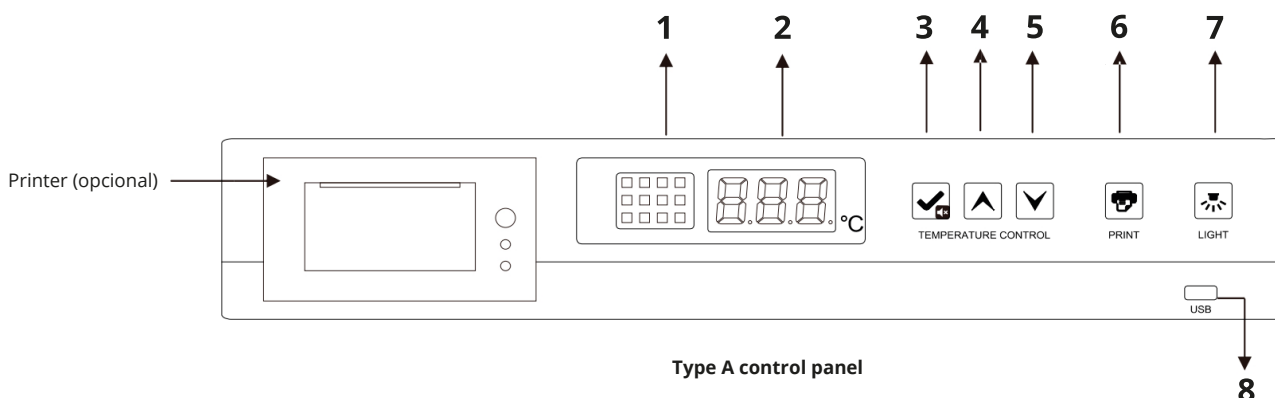

1. Function description od Type A control panel (except TH-081)

1) Description of display icon indicator (as shown in the right diagram);

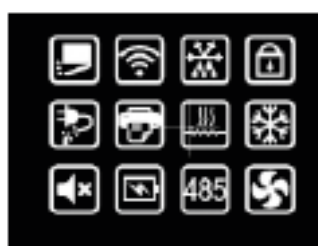Door switch	WIFI	Defrosting	Key lock
Power failure	Print	Door heating	Refrigeration
Mute	Electric quantity	Serial port	Fan

a Door switch indicator

When the door is opened, the door switch indicator is on; when the door is closed, the door switch indicator is off.

b WIFI indicator

WIFI pilot lamp is on when WIFI is connected; it goes out when WIFI is disconnected.

c Defrosting indicator

When the refrigerator enters the defrosting state, the defrosting indicator is always on; When the refrigerator exits the defrosting state, the defrosting indicator is always off.

d Key lock indicator

When the keys are in the lock state, none of the keys respond and the key lock lights up. In this case, after pressing the up key+down key for 3s, you will be prompted to enter the password, which is defaulted as "005". After entering it correctly, press the Set/Mute Multiplex key and the key lock will be released. At this time, the key lock indicator will be off. In the unlocked state, if no key is pressed for 60 seconds, the key lock is started and the key lock indicator is on. Long press the key + down key for 3 seconds, and the keys are locked.

e Power-ff indicator

The refrigerator is normally powered by 220V/110V. When the input power is turned off, the buzzer will be triggered, the digital tube flashes the power off code "PF" alternately at 3s intervals and the power-off indicator is on. When the input power is turned on, it returns to normal and the power-off indicator is off.

f Print indicator

When the printer is not working, the indicator is off; When the printer is working, the indicator is on.

g Door heating indicator

When the door heating is turned off, the door heating indicator is off; When the door heating is turned on, the door heating indicator is on.

h Refrigeration indicator

If the compressor is in working condition, the refrigeration indicator is on; if the compressor is in shutdown state, the colling indicator is off.

i Mute indicator

When the alarm tone key is muted, the indicator is on; When the alarm mute function is cancelled, the indicator is off.

j Low battery level indicator

When the battery voltage is less than 8V, the buzzer will be triggered, the low battery indicator will be on and the digital tube will flash the low battery code "BL" alternately at 3s intervals; When the battery voltage is greater than 12V, the buzzer will be turned off, the low battery indicator will be off and the digital tube will resume normal display.

k Serial port indicator

When the reserved RS-485 serial port is not connected to the equipment, the serial port indicator will be off; When the equipment is successfully connected to the reserved RS-485 serial port, the serial port indicator will be on.

l Fan indicator

When the evaporator fan is turned on, the fan indicator is on; When the evaporator fan is turned off, the fan indicator is off.

2) Temperature display window, which displays the average temperature inside the cabinet in °C under normal operating;

View ambient temperature:

In the key lock state, press  the key and digital tube displays the ambient temperature and returns to normal display after 5s without key operation or pressing  and . In the key unlocked state, press  and digital tube displays ambient temperature and returns to normal display after 5 seconds without key operation.

Humidity check: key unlocked state, long press and , humidity of the digital tube display, do not press any key operation after 5 seconds or press  and  to return to normal display.

3)  is the set/mute key;

In case of no alarm state and key unlock status, press  and display the ambient temperature for 5s and then return to normal display; In unlock status, press  for more than 3 seconds and enter the user menu.

When the buzzer is triggered (including cabinet high temperature alarm, door opening alarm, sensor failure alarm, etc.) and in the key unlock state, press  for the first time, and the buzzer stops ringing, and the ambient temperature is displayed for 5 seconds, after which the normal display resumes (pressing the mute button is only to turn off the buzzer for alarming this abnormal state, for example, troubleshooting, and the buzzer will be triggered next time for any abnormality). Then press  again, trigger the buzzer, display the ambient temperature for 5 seconds, and resume the display of the chamber temperature and alarm state. In the key unlocked state,  can be used as a configuration key.

In the unlock state and parameter setting mode, press this key to display the values and parameter names. If the pressing time is longer than 3 seconds, save the settings and return to the normal interface.

4)  is a up key;

In parameter setting mode, move to the next parameter or increase the parameter value. For example, when setting the set temperature, increase the set temperature value. When setting the parameter value, press the up button and the parameter will increase rapidly. Under normal conditions, long press the up for 3 seconds to import the data of the USB flash drive in 12 months.

5)  is a down key;

In parameter setting mode, move to the previous parameter or decrease the parameter value. For example, when setting the set temperature, reduce the set temperature. When setting the parameter value, long press the down button and the parameter will decrease rapidly.

6)  is a print key;

The system can keep 7 days of data for printing and press the print key to print the temperature within the set time period.

7)  is a light switch key

When the equipment is powered on, the lights are turned off by default.

8) USB data export;

Automatic export: When the U disk is connected to the USB interface, the recorder buzzer will chirp once and display "On". PDF files of data that not currently exported will be generated in the U disk. After data transmission, the buzzer will chirp once more again and display "End". After 6s, it will return to normal display.

Note: When there is less data, the digital tube will not display "on" and "End".

Manual export: In the key unlock state, and when the USB flash drive is connected and the file is not being generated, press the key up for 3 seconds and the digital tube will display "d01". Press the up key or down key to set "d00 ~ d12" and press  key to obtain the file generation (d00) or generate the PDF file of the record data of the previous months (1-12).

2. Function setting of Type A control panel

1) After powering on, the equipment can enter the working state;


2) User parameter setting:

Unlock: Under normal operating state, simultaneously press  and  keys for 3s and the digital tube will display the parameter code "0000"; Pressing to enter the password "0005" and holding to unlock.

Then press  the key for 3s, the digital tube will display the parameter code "PS1" and enter the setting and adjustment parameters.

a Use the  or  key to scroll the parameters;

b Press  the key to display the corresponding parameter value;

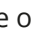c Use the  or  key to scroll through the parameters;

d Use  to  temporarily store the modified values and return to the display parameters;

e If other parameters are modified, repeat steps 1 to 4

f Press  for more than 3 s, save the modified parameters and return to the display parameters category.

3) Press  for more than 3 s, or press no key within 60 S to exit the parameter setting program.

4) Parameter display

No.	Menu item	Parameter Range	Suggested settings	Remarks
1	MAX.	-	-	The highest temperature since last clearance
2	MIN.	-	-	The lowest temperature since last clearance
2	CLR	-	-	Clearance of the Max and Mini temperature records
4	Set	0.0 - 10.0	5.0	Temperature setting
5	H	0.0 - 10.0	5.0	Set value of high temperature alarma set+H;When H=0, High temp alarm is disabled; When the alarm is over high temp set, H1 will be displayed on the controller
6	L	0.0 - 10.0	5.0	Set value of low temperature alarma set-L;When L=0, Low temp alarm is disabled; When the alarm is below low temp set, L1 will be displayed on the controller
7	n	Set logger module time - year	-	-
8	y	Set logger module time - month	-	-
9	r	Set logger module time - day	-	-
10	s	Set logger module time - hour	-	-
11	F	Set logger module time - minute	-	-
12	Pt	0-240 min	20	-
13	tH1	20.0-50.0 °C	40.0	-

No.	Menu item	Parameter Range	Suggested settings	Remarks
14	P1	1. Automatic heating mode 1 2. Automatic heating mode 2 3. Automatic heating mode 3 4. Always on 5. Always off	1. (Set to 4 when the door is with condensation) (Set to 1 when the door is a glass door) (Set to 5 when the door is a foaming door)	Mode 1: It shall be judged as once after the door is opened and closed once and heated for 5min (time setting). If the door is opened and closed again during the heating period, the heating time will be updated again; Mode 2: When the compressor runs, heater is on; When the compressor stops, the heater will automatically off after one minute. Mode 3: When the humidity in the cabinet is more than 80%, the door heating is on, and when the humidity in the cabinet is moderately less than 60%, the door heating is off; Mode 4: Door heating is always on; Modo 5: Door heating is always off.
15	Ps1	0000-9999	0005	User menu password settings
16	b1	-	-	Repair Information 1
17	b2	-	-	Repair Information 2

Quick setting of time after power on


After the power-on self-test on the display board is completed, the quick setting menu is displayed

Menu item	Menu	Menu description	Set range	Default	Unit
Quick Setting Menu	n	Set logger time - year	10-50		/
	y	Set logger time - month	1-12	--	/
	r	Set logger time - day	01-31	--	/
	S	Set logger time - hour	00-23	--	/
	F	Set logger time - minute	00-59	--	/
	Pt	Print interval	0-240	20	min
	SCY	Temperature data recording period	0-240 0: Shielded recorder	10	min

If there is no operation for 60 seconds under the quick setting menu, it will automatically exit the quick setting menu and return to normal display.

3. Alarm display

Code	Error Description
H1	High temperature alarm
L1	Low temperature alarm
H2	Alarm for high ambient temperature
H3	Condenser overheat alarm
do	Door opening alarm
PF	Power failure alarm
bL	Battery low alarm
Er	The recorder is not connected.
LoF	The recorder did not start
EE.	Communication failure

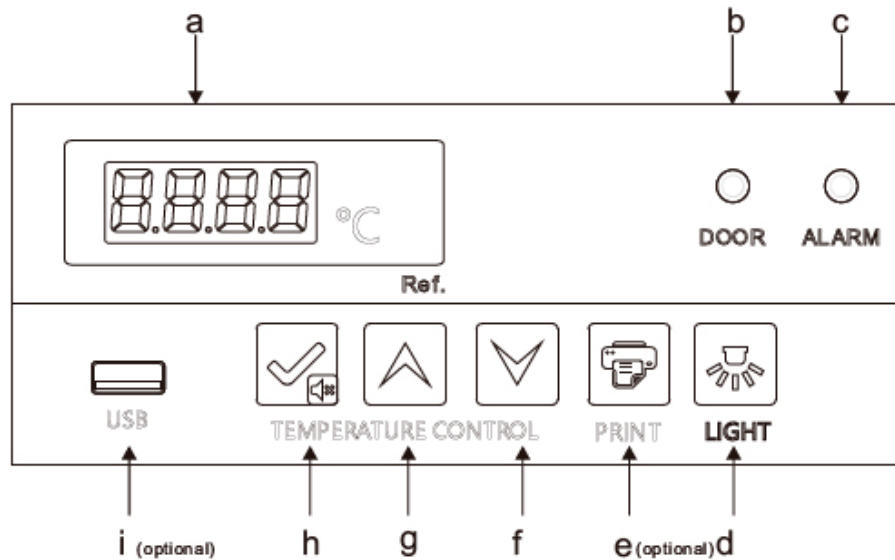1. Function description of Type B control panel (Applicable to TH-081)

- a** It is temperature display window, which displays the average temperature inside the cabinet in °C in under normal operation; different prompt characters can be displayed in the setting state (see below for details).
- b** Door open indicator: When the refrigerator door is opened, the indicator lights up. After more than 1 minute, the door opening alarm indicator will be on and the buzzer will triggered, displaying "do".
- c** Fault indicator: When the product runs normally, the indicator is off; The indicator is on for operation abnormality.
- d** "LIGHT": After the machine is powered on, the light is off by default and the on/off of the light can be adjusted with the on/off key.
- e** "PRINT" (optional): The system can keep 7 day of data for printing. Press the print key to print the temperature within the set time period.
- f** : In parameter setting mode, reduce the parameter value. For example, when setting the temperature, reduce the set temperature. When setting the parameter value, long press the down button and the parameter will decrease rapidly.
- g** : In parameter setting mode, increase the parameter value. For example, when setting the temperature, increase the set temperature value. When setting the parameter value, long press the up button and the parameter will increase rapidly. Under normal conditions, long press the up key for 3 seconds to import the data from the USB flash drive in 12 months.
- h** is set/silence key; In case of no alarm state and key unlock state, press , and display the ambient temperature for 5 seconds and then return to normal display; In unlock status, press for more than 3s and enter the user menu.
When buzzer is triggered (including cabinet high alarm, door opening alarm, sensor failure alarm, etc.) and button unlock state of the, press for the first time, and the buzzer stops ringing, and the ambient temperature is displayed for 5 seconds, after which normal display is resumed (pressing the mute button is only to turn off the buzzer to alarming this abnormal state, for example, trouble removal, and the buzzer will be triggered next time for any abnormality). Then press again, trigger the buzzer, display the ambient temperature for 5 seconds, and then resume the display of cabinet temperature and alarm state. In the key unlocked state, can be used as a setting key.
In the unlock state and parameter setting mode, press this key to display the values and parameter names. If the pressing time is longer than 3 seconds, save the setting and return to normal interface.
- i** USB interface. Automatic USB data export: When the USB interface is connected to the USB flash drive, the buzzer of the recorder beeps once, displaying "ON", and PDF files of the data of the current month and previous month are generated in the USB flash drive. After the data transmission is completed, the buzzer beeps once, displaying "End", and the normal display resumes after 6 seconds.

⚠ Note: When there is less data, "on" and "end" prompts are not displayed

Manual export of USB data: In the key unlock state, when the USB flash drive is connected and the file is not being generated, press the up key for 3 seconds and the digital tube in the lower chamber will display "d01". Press the key up or down key to set "d00~d12" and press key to cancel the file generation (d00) or generate the PDF file of the record data of the previous month (1-12).

Note: When the alarm of the digital tube flashes and displays "LoF", the recorder is not started; Meanwhile press **↔** and **↕** keys for 3s, and "LoF" disappears, the recorder starts.

2. Function setting of Type B control panel

1) After powering on, the equipment can enter the working state;

2) User parameter settings:

Unlock: normal operating state, simultaneously press **↔** and **↕** keys for 3s and the digital tube will display the parameter code "0000"; By pressing **↔** to enter the password "0005" and holding **↕** to unlock. Then press **↕** key for 3s, the digital tube will display the parameter code "PS1" and enter the setting and adjustment parameters.

- a Use **↔** and **↕** key to scroll the parameters;
- b Press **↕** key to display the value of the corresponding parameter;
- c Use **↔** or **↕** key to scroll through the parameters;
- d Use to temporarily store the modified values and return to the display parameters;
- e If other parameters are changed, repeat steps 1 - 4.

3) Press **↕** for more than 3 s, save the modified parameters and return to the display parameters category.

4) Press **↕** for more than 3 s, or press no key within 60 S to exit the parameter setting program.

3. Parameter display

No.	Menu item	Parameter Range	Suggested settings	Remarks
1	MAX.	-	-	The highest temperature since last clearance
2	MIN.	-	-	The lowest temperature since last clearance
2	CLR	-	-	Clearance of the Max and Mini temperature records
4	Set	0.0 - 10.0	5.0	Temperature setting
5	H	0.0 - 10.0	5.0	Set value of high temperature alarm set+H;When H=0, High temp alarm is disabled; When the alarm is over high temp set, H1 will be displayed on the controller
6	L	0.0 - 10.0	5.0	Set value of low temperature alarm set-L;When L=0, Low temp alarm is disabled; When the alarm is below low temp set, L1 will be displayed on the controller
7	n	Set logger module time - year	-	-
8	y	Set logger module time - month	-	-
9	r	Set logger module time - day	-	-
10	s	Set logger module time - hour	-	-
11	F	Set logger module time - minute	-	-
12	Pt	0-240 min	20	-
13	tH1	20.0-50.0 °C	40.0	-

No.	Menu item	Parameter Range	Suggested settings	Remarks
14	P1	1. Automatic heating mode 1 2. Automatic heating mode 2 3. Automatic heating mode 3 4. Always on 5. Always off	1. (Set to 4 when the door is with condensation) (Set to 1 when the door is a glass door) (Set to 5 when the door is a foaming door)	Mode 1: It shall be judged as once after the door is opened and closed once and heated for 5min (time setting). If the door is opened and closed again during the heating period, the heating time will be updated again; Mode 2: When the compressor runs, heater is on; When the compressor stops, the heater will automatically off after one minute. Mode 3: When the humidity in the cabinet is more than 80%, the door heating is on, and when the humidity in the cabinet is moderately less than 60%, the door heating is off; Mode 4: Door heating is always on; Modo 5: Door heating is always off.
15	Ps1	0000-9999	0005	User menu password settings
16	b1	-	-	Repair Information 1
17	b2	-	-	Repair Information 2

Quick setting of time after power on

After the power-on self-test on the display board is completed, the quick setting menu is displayed

Menu item	Menu	Menu description	Set range	Default	Unit
Quick Setting Menu	n	Set logger time - year	10-50		/
	y	Set logger time - month	1-12	--	/
	r	Set logger time - day	01-31	--	/
	S	Set logger time - hour	00-23	--	/
	F	Set logger time - minute	00-59	--	/
	Pt	Print interval	0-240	20	min
	SCY	Temperature data recording period	0-240 0: Shielded recorder	10	min

If there is no operation for 60 seconds under the quick setting menu, it will automatically exit the quick setting menu and return to normal display.

Code	Error Description
H1	High temperature alarm
L1	Low temperature alarm
H2	Alarm for high ambient temperature
H3	Condenser overheat alarm
do	Door opening alarm
PF	Power failure alarm
bL	Battery low alarm
Er	The recorder is not connected.
LoF	The recorder did not start
EE.	Communication failure

6.2 Optional function

6.2.1 Printer


The printer has been installed with a paper roll at the factory. If the paper roll runs out after a long period of use, you can buy the same paper roll (size: thermal paper, paper width: 57.5 ± 0.5 mm, the outer diameter of the reel: no more than 40 mm, i.e., rear roll paper RM57 * 40).

Description of printer panel:

Open button. Press to open the cover

SEL button, indicator light, for factory setting. Do not press it;

Lf button. The green indicator is the power indicator and normally lit when the power is turned on;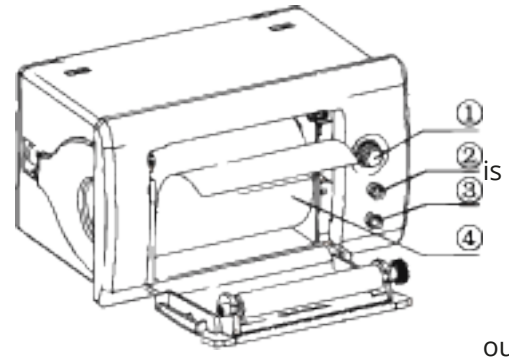


Install the paper roll:

Press the open button 1, open the cover, install the paper roll, close the cover plate and allow the head of the roll paper to slightly extend of the cover plate.

6.2.2 Remote alarm terminal, RS485 interface and network interface


Remote alarm terminal, RS485 interface and network interface are installed in the lower part of the back of the freezer body.

Note: These configurations are standard for some models.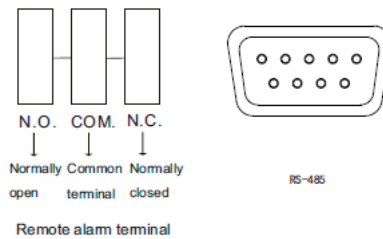

6.2.3 Chart recorder

This is an optional function. If the function is selected, refer to the "Chart recorder operating instructions" in the accessory bag.

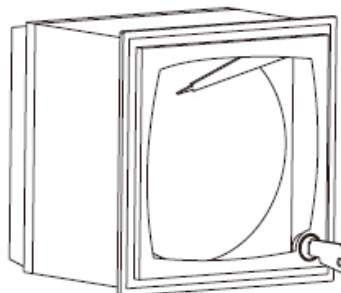7. Maintenance and Service

Warning:

- In order to prevent people from getting electric shock or injury, please cut off the power supply of the equipment before any repairs and maintenance.

- Make sure you don't inhale drugs or suspended particles around you when servicing the equipment, otherwise it will damage your health.

7.1 Equipment Maintenance

Defrosting: The equipment will automatically defrost during the working process for convenience of use.

Cleaning and maintenance: The equipment should be cleaned and maintained regularly (for safety, please unplug the power plug), and the inner and outer surfaces of the cabinet should be wiped with a warm, damp soft cloth. Condenser cleaning: clean the dust on the condenser regularly to avoid affecting the cooling effect (cleaning method: first open the screws on the ventilation hood or front hood and use a brush to gently brush away the dust on the surface of the condenser).

Notes:

- Do not sprinkle water directly on the cabinet, lest the insulation performance of electrical components decrease and metal parts rust.
- Do not use hot water, corrosive detergent or organic solvent to clean the cabinet.
- Do not place heavy objects on top of the equipment, as the equipment may deform under pressure.
- The integrated overload protector within our products is programmed to shut down the power supply automatically within respect short circuits and/or system overload for the sake of protecting the entire system and crucial components such as the compressor.

7.2 Equipment Discontinuation

Deactivation: If the equipment is stored in an unsupervised area for a long time and not used, the power supply should be cut off and the inner and outer surfaces of the cabinet should be cleaned with warm, damp soft cloth aired and sealed. The equipment must be locked to ensure that children cannot open the cabinet door.

Scrapping: When the equipment reaches the end of its service life, it should be scrapped and must be handed over to a qualified professional recycling agency for disposal per local regulations. Non-professionals are not allowed to disassemble or break down the equipment without authorization. The scrapped equipment should be placed in a designated area inaccessible to children to avoid danger.

7.3 Maintenance, Replacement and Recovery of Rechargeable Batteries

Battery installation position: top of the cabinet and bottom inside the electrical box.

1. Battery maintenance

- 1 In order to prolong the service life and avoid the product being left unused for a long time, the product must be operated for more than 24 hours every month to facilitate charging.
- 2 The equipment should be connected to the power supply regularly (generally once a month) and turn on the power switch and battery switch for a period of time to complete charging if it is not used for a long time, and each charging time should not be less than 24 hours;
- 3 Power switch and battery switch should be turned off in time after discontinuation of equipment, otherwise it will cause battery capacity loss, and in serious cases, battery permanent damage could happen;
- 4 The main power switch must be turned off during long-term power outage or during transportation; otherwise, long-term discharge causes power loss in the battery and even permanent damage to the battery, and the display is abnormal after re-energizing.
- 5 The battery is a consumable, with a service life of about 2-3 years. If the battery is improperly, such as power loss or reaching the end of battery life, a low battery alarm (refrigeration is not affected, but there is an alarm failure and influence to the use function of the printer, please contact the after-sales staff of local distributors for replacement).

2. Battery replacement and recovery

- 1 Turn off the power switch and pull out the power plug from the outlet;
- 2 Remove the screws on the electrical box with a screwdriver. (Note: There are high-voltage electrical components in the electrical box, therefore, before opening, turn off the power supply, unplug the power plug and turn off the power lock switch of the refrigerator, and the electrical box can only be opened by

qualified engineers or maintenance personnel);

- 3 Pull out the battery connecting cable; (Before unplugging the battery fixing cable, pay attention to the cathode/anode of the battery and the sequence of connection cables, so as to prevent burning down the control system due to the installation of the new battery with cathode/anode reversed. Generally, the red wire is connected to the cathode and the black wire is connected to the anode);
- 4 Use a screwdriver to remove the fastening screws on the battery fixing plate and take out the battery;
- 5 For a replaced battery that can be recycled, please contact the local battery recycling agency.

Tips:

In order to effectively ensure that the battery replacement meets the requirements of the control system and avoid the influence of improper operation of the system, please contact our after-sales personnel for replacement or guidance.

When the circuit is short circuit or overloaded, the overload protector will disconnect the power supply of the device. Please contact after sales at this time.

8. Troubleshooting and Maintenance Services

Any product may fail. Please observe the operation of equipment in time during use. If there is any abnormality, please check and handle it according to the following table. If the abnormality can't be changed, please inform our service center in time, and we will serve you wholeheartedly to avoid losses.

Problems	Causes and solutions
The equipment does not work	Make sure the power outlet is energized. Make sure that the power plug is plugged in and not loose. Make sure that the power fuse is not disconnected. Make sure that the supply voltage is adequate, neither too low nor too high.
Compressor is not running	Please make sure that the temperature is set correctly. Check if the temperature inside the cabinet is too low.
Temperature does not reach the set value	Make sure the door is tightly closed and do not open it too many times in a short period of time. Please do not put too many items at a time. Make sure that the ambient temperature is not too high.
High noise	Make sure the cabinet is placed on a flat surface. Make sure that the cabinet does not make contact with the wall.
Condensation on the cabinet surface	In rainy and humid seasons, condensation on the door is normal and should be wiped off with a dry cloth.
Undesirable odor	Equipment needs to be cleaned. Items with strong odor are not packed.
The door is not properly closed and cold air leaks in.	Please make sure that the temperature is set correctly. Check if the temperature inside the cabinet is too low.
The alarm lamp flashes and the buzzer is activated.	After using the equipment for a period of time, the door seal becomes hard and deformed. Maintenance method: Blow the deformed part of the hot seal with a blower to soften it, and then close and compress it after the door seal softens.
Lighting damage (refrigeration equipment)	Call the after-sales hotline and contact Thermolabil® after-sales service personnel for replacement. Do not replace parts yourself.

The following conditions are not faults:

- 1 When the compressor starts and stops, the equipment parts will make a slight impact sound;
- 2 When overheated items are put in after the door is opened, high temperature alarm and high humidity alarm will be triggered in the control system (if this function is available, please refer to the alarm display table);

Solution: Place items into the equipment after they are naturally cooled. Put the items in small quantities in batches and do not put too many at a time. After the system runs stably, the high temperature alarm and high humidity alarm will be released automatically.

3 Slight sound of running water caused by refrigerant flow in the pipeline.

 **Warning**

- The system contains refrigerant under high pressure. Do not tamper with the system. It must be serviced by suitably qualified persons only.
- Connect to the potable water supply only.
- Do not store explosive substances such as aerosol cans with a flammable propellant in this appliance.
- If the symbol ISO 7000 1701 (2004-01) is used, its meaning will be explained.
- In order to reduce flammability hazards the installation of this appliance must only be carried out by a suitably qualified person.
- The refrigeration system is under high pressure. Do not tamper with it. Contact qualified service personnel before disposal.

 **Notes:**

- The equipment can only be repaired, maintained or improved by engineers certified by Thermolabil®, so as to ensure the normal operation of the equipment and compliance with the corresponding safety standards.
- Please clean and disinfect the equipment before notifying the maintenance engineer; During the warranty period of the equipment, the Company will not undertake the warranty obligation if the fault or damage is caused by improper use of the user.
- Ambient temperature for storage: -40°C +55°C, relative humidity: 10% ~ 90%.

9. Specifications

Model	Ambient temperature (°C)	Climate type	Refrigerant and loading amount	(Rated) voltage (V~)	Rated frequency (Hz)	Temperature range (°C)	Volume (L)	(Rated) current (A)	Weight (kg)	Exterior dimensions (DxWxH) (mm)
TH-076	16~32	N	R600a/18g	110	60	2~8	76	1,24	41	566x542x764
TH-081	16~32	N	R290/26g	110	60	2~8	130	2,28	51	625x650x810
TH-176	16~32	N	R290/26g	110	60	2~8	315	3,83	87	673x650x1762

10. Packaging List

Item	Operation Manual	Certificate of conformity	Accessory package	Key
Quantity		1	1	2

*The specific packing list shall be subject to the physical objects received.